

Application Note

Downloading a Method to the i-LAB®

You have received a customized method from micrOptix Technical Support by e-mail and need to transfer this method to your i-LAB®. These steps will outline how to import the method into the Spectrum Software and download it into the i-LAB®.

- 1) Save the file to disk or desktop by dragging the .xml attachment from the e-mail to your desktop.
- 2) Connect the i-LAB® with the USB cable and power on the i-LAB®.
- 3) Open the **i-LAB® Spectrum** software and from the Menu in the left bar select to open **Method Designer**.
- 4) When open, select the **Sharing** tab, then select **Import from File**.
- 5) A window will open, select **Desktop** then select the **.xml file** you just saved to your desktop.
- 6) Select **Open**.
- 7) Click on the "Send to i-LAB®" tab.
- 8) File will download to your i-LAB® **Methods** folder.
- 9) When finished, detach the i-LAB® USB cable.
- 10) Consult the method instructions for details on running the method.

Thank You for contacting micrOptix Technical Support.

micrOptix Technologies, LLC

284 Main Street
Section L Suite 400
Wilton Maine 04294
Tel: 207-645-3600
Fax: 207-645-3633
Email: TechSupport@microSpectralSensors.com

