

i-LAB Spectrum: Windows Vista Software Installation Bulletin

Issue: i-LAB Spectrum program may give an error message stating the “*database is read-only*” when installed in a Windows Vista environment.

Cause: This is caused by the enhanced security settings in Vista, causing permission violations on files and folder.

Fix: To fix the problem, perform the following:

- 1) Navigate to the iLAB/Database folder. By default this is installed in C:\Program Files\MicroSpectral Sensors.
- 2) Right-click on the iLab/**Database** folder
- 3) Click **Properties**
- 4) In the Database properties window, click the **Security** tab
- 5) Under the Group or User Names, click the **Edit** button.
- 6) Select the appropriate user that is logged in, or the “**Users**” group in the grid
- 7) Under Permissions for User, Check “**Allow**” under **Full Control**, and under **Modify**
- 8) **Re-boot the PC.**

